# Product Catalogue 2011 - 2012

**DNA/RNA Purification Kits**

## GeneON GmbH

Hubertusstraße 20  
D-67065 Ludwigshafen

Phone: +49-621-5720 864

Fax: +49-621-5724 462

E-Mail: [info@geneon.net](mailto:info@geneon.net)

Web: [www.geneon.net](http://www.geneon.net)

## Content

### DNA Purification Kits

Bacterial DNA Extraction Kit (Proteinase K included) .....	8- 9
Blood DNA Extraction Kit (Proteinase K included) .....	10-11
Plant DNA Extraction Kit (Proteinase K included) .....	12-13
Tissue DNA Extraction Kit (Proteinase K included) .....	14-15
Plasmid DNA Extraction Kit (Proteinase K included).....	16-17
Food DNA Extraction Kit (Proteinase K included) .....	18-19
PCR Clean-up Kit.....	20-21
Gel DNA Recovery Kit.....	22-23
AmbiClean Kit (PCR & Gel).....	24-25

### RNA Purification Kits

Total RNA Extraction Kit (Proteinase K & DNase I included) .....	28-29
Blood Total RNA Extraction Kit (Proteinase K & DNase I included) .....	30-31
Viral Nucleic Acid Extraction Kit (Proteinase K, DNase I & Carrier RNA included) .....	32-33

# DNA/RNA Purification Kits

DNA Purification Kits ..... 7

RNA Purification Kits .....27

## DNA Purification Kits

- Bacterial DNA Extraction Kit (Proteinase K included) ..... 8- 9
- Blood DNA Extraction Kit (Proteinase K included) ..... 10-11
- Plant DNA Extraction Kit (Proteinase K included) ..... 12-13
- Tissue DNA Extraction Kit (Proteinase K included)..... 14-15
- Plasmid DNA Extraction Kit (Proteinase K included)..... 16-17
- Food DNA Extraction Kit (Proteinase K included) ..... 18-19
- PCR Clean-up Kit ..... 20-21
- Gel DNA Recovery Kit ..... 22-23
- AmbiClean Kit (PCR & Gel) ..... 24-25

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- Columns
- Collection tubes
- Resuspension Puffer 1
- Resuspension Puffer 2
- Bacterial Genomic Binding Puffer
- Waschpuffer
- Elutionspuffer
- Proteinase K Lösung \*
- Handbuch

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# Bacterial DNA purification

## Description:

The GF-1 Bacterial DNA Extraction Kit is designed for rapid and efficient purification up to 20µg of a high molecular weight genomic DNA from 1-3 ml either Gram-negative or Gram-positive bacteria. This kit uses a specially-treated glass filter membrane fixed into a column to efficiently bind DNA in the presence of high salt.

This bacterial purification kit applies the principle of a minicolumn spin technology and the use of optimized buffers to ensure that only DNA is isolated while cellular proteins, metabolites, salts and other low molecular weight impurities are removed during the subsequent washing steps.

High-purity genomic DNA is eluted in water or low salt buffers and has a A260/280 ratio between 1.7 and 1.9 making it ready to use in many routine molecular biology applications such as restriction enzyme digestion, Southern blotting, PCR, DNA fingerprinting and other manipulations.

## Kit components:

- Columns
- Collection tubes
- Resuspension Buffer 1
- Resuspension Buffer 2
- Bacterial Genomic Binding Buffer
- Wash Buffer
- Elution Buffer
- Proteinase K solution\*
- Handbook

\* Please refer to Reconstitution of Solutions and Storage and Stability before using this kit.

Cat.-no	Description	Amount
BA05	Bacterial DNA extraction kit	5 preps
BA050	Bacterial DNA extraction kit	50 preps
BA100	Bacterial DNA extraction kit	100 preps

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# Blood DNA Purification

## Description:

The GF-1 Blood DNA Extraction Kit is designed for rapid and efficient purification of genomic DNA from fresh anti-coagulated whole blood and frozen white blood cell. The purification is based on the usage of denaturing agents to provide lysis of cells, denaturation of proteins and subsequently release of genomic DNA. Special buffers provided in the kit are optimized to enhance binding of DNA onto a specially-treated glass filter membrane for efficient recovery of highly pure genomic DNA.

## Features:

- Purification of up to 1mL of whole blood
- Yields up to 20µg of DNA
- Mini-column spin technology
- No organic-based extraction required
- Highly pure genomic DNA ready to use for routine molecular biology applications such as restriction enzyme digestion, PCR, Southern blotting, DNA fingerprinting, etc.

## Application:

For high quality blood genomic DNA isolation

## Kit components:

- GF-1 columns
- Collection tubes
- Blood Lysis Buffer (Buffer BB)
- Wash Buffer 1 (concentrate)\*
- Wash Buffer 2 (concentrate)\*
- Elution Buffer
- Proteinase K\*
- Handbook

\* Please refer to Reconstitution of Solutions and Storage and Stability.

Cat.-no	Description	Amount
BD05	Blood DNA extraction kit	5 preps
BD050	Blood DNA extraction kit	50 preps
BD100	Blood DNA extraction kit	100 preps

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- GF-1 columns
- Collection tubes
- Plant Tissue Lyse Puffer (Puffer PL)
- Plant Genomic Binding Puffer (Puffer PB)
- Waschpuffer (Konzentrat)\*
- Elutionspuffer
- Proteinase K\*
- Handbuch

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# Plant DNA Purification

## Description:

The GF-1 Plant DNA Extraction Kit is designed for rapid and efficient purification of genomic DNA from a variety of plant tissues without the need for precipitation or organic extraction. This kit uses a specially treated glass filter membrane fixed into a column to efficiently bind DNA in the presence of high salt. The kit applies the principle of a minicolumn spin technology and the use of optimized buffers to ensure that only DNA is isolated while proteins and other impurities, additives, preservatives are removed during the subsequent washing steps. High purity genomic DNA is then eluted in water or low salt buffers has a A260/280 ratio between 1.7 and 1.9, making it ready to use in many routine molecular biology applications such as restriction enzyme digestion, Southern blotting, DNA fingerprinting, PCR and other manipulations

## Features:

- Yields up to 20µg of DNA
- Mini-column spin technology
- No organic-based extraction required

## Application:

For high quality plant genomic DNA isolation

## Kit Components:

- GF-1 columns
- Collection tubes
- Plant Tissue Lysis Buffer (Buffer PL)
- Plant Genomic Binding Buffer (Buffer PB)
- Wash Buffer (concentrate)\*
- Elution Buffer
- Proteinase K\*
- Handbook

\* Please refer to Reconstitution of Solutions and Storage and Stability.

Cat.-no	Description	Amount
PT05	Plant DNA extraction kit	5 preps
PT050	Plant DNA extraction kit	50 preps
PT100	Plant DNA extraction kit	100 preps

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- GF-1 columns
- Collection tubes
- Tissue Lyse Puffer (Puffer TL)
- Lyse Verstärker
- Tissue Genomic DNA Binding Puffer (Puffer TB)
- Waschpuffer (Konzentrat)\*
- Elutionspuffer
- Proteinase K\*
- Handbuch

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# Tissue DNA Purification

## Description:

The GF-1 Tissue DNA Extraction Kit is designed for rapid and efficient purification of genomic DNA from various tissue samples such as kidney, heart, lungs, brain, muscles, liver, spleen, etc. The purification is based on the usage of denaturing agents to provide lysis of tissue cells, denaturation of proteins and subsequently release of genomic DNA. Special buffers provided in the kit are optimized to enhance the binding of DNA onto a specially-treated glass filter membrane for efficient recovery of highly pure genomic DNA.

## Features:

- Yields up to 20µg of DNA
- Mini-column spin technology
- No organic-based extraction required
- Highly pure genomic DNA ready to use for routine molecular biology applications such as restriction enzyme digestion, PCR, - Southern blotting, DNA fingerprinting, etc.

## Application:

For high quality tissue genomic DNA isolation

## Kit Components:

- GF-1 columns
- Collection tubes
- Tissue Lysis Buffer (Buffer TL)
- Lysis Enhancer
- Tissue Genomic DNA Binding Buffer (Buffer TB)
- Wash Buffer (concentrate)\*
- Elution Buffer
- Proteinase K\*
- Handbook

\* Please refer to Reconstitution of Solutions and Storage and Stability.

Cat.-no	Description	Amount
TD05	Tissue DNA extraction kit	5 preps
TD050	Tissue DNA extraction kit	50 preps
TD100	Tissue DNA extraction kit	100 preps

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# Plasmid DNA Purification

## Description:

The GF-1 Plasmid DNA Extraction Kit is designed for rapid and efficient purification of high copy and low copy plasmid DNA without the need for precipitation or organic extractions. This it uses a specially-treated glass filter membrane fixed into a column to efficiently bind DNA in the presence of high salt. Combining alkaline lysis-SDS and minicolumn spin technology, up o 20µg of plasmid DNA from bacterial cultures can be isolated. Multiple samples can be processed rapidly and with practice, the purification takes less than 30 minutes. Optimized buffers ensure only highly pure plasmid DNA is extracted and is ready to use in many routine molecular biology applications such as restriction enzyme digestion, radioactive/fluorescence DNA sequencing, PCR, ligation, transformation and other manipulations.

## Features:

- Yields up to 20µg of DNA
- Multiple samples can be processed rapidly in less than 30 minutes
- No organic-based extraction required

## Application:

For high quality plasmid DNA isolation

## Kit Components:

- GF-1 columns
- Collection tubes
- Solution 1 (S1)\*
- Solution 2 (S2)\*
- Neutralizing Buffer (Buffer NB)
- Wash Buffer (concentrate)\*
- RNase A (DNase-free)\*
- Elution Buffer
- Handbook

\* Please refer to Reconstitution of Solutions and Storage and Stability before using this kit.

Cat.-no	Description	Amount
PL05	Plasmid DNA extraction kit	5 preps
PL050	Plasmid DNA extraction kit	50 preps
PL100	Plasmid DNA extraction kit	100 preps
PL200	Plasmid DNA extraction kit	200 preps

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## Description:

The GF-1 Food DNA Extraction Kit is designed for rapid and efficient purification of genomic DNA from up to 100mg of raw or processed food from plant, animal or mixed origins. This kit uses a specially-treated glass filter membrane fixed into a column to efficiently bind DNA in the presence of high salt. The kit applies the principle of a minicolumn spin technology and the use of optimized buffers to ensure that only DNA is isolated while proteins and other impurities, additives, preservatives are removed during the subsequent washing steps. High purity genomic DNA is then eluted in water or low salt buffers, ready to use in many routine molecular biology applications such as PCR and restriction enzyme digestion. This kit is suitable for use in GMO testing and has been tested on various raw samples such as soybean, maize and processed foods (soy flour, soymilk, soy sauce, sausage, meat, cereal meal, and cocoa).

## Features:

- Mini-column spin technology
- No organic-based extraction required
- Highly pure DNA, ready to use for routine molecular biology applications such as PCR

## Application:

For purification of DNA from raw or processed food

## Kit Components:

- GF-1 columns
  - Collection tubes
  - Food Lysis Buffer (Buffer FL)
  - Food DNA Binding Buffer (Buffer FB)
  - Wash Buffer 1 (concentrate)\*
  - Wash Buffer 2 (concentrate)\*
  - Elution Buffer
  - Proteinase K\*
  - Handbook
- \*Please refer to Reconstitution of Solutions and Storage and Stability before using this kit.

Cat.-no	Description	Amount
TD05	Food DNA extraction kit	5 preps
TD050	Food DNA extraction kit	25 preps

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## PCR Clean-up Purification

### Description:

The GF-1 PCR Clean-up Kit is designed for rapid and efficient clean up of DNA ranging from 100bp to 20kb. The Kit contains special buffers to provide the correct salt concentration and pH for efficient recovery (80 - 90%) of DNA. This kit uses a specially treated glass filter membrane fixed into a column to efficiently bind DNA in the presence of high salt. It applies the principle of a mini-column spin technology and is well suited for the removal of excess dNTPs, short oligo fragments, mineral oil, enzymes from a PCR reaction product, proteins after restriction enzyme treatment and dephosphorylation and residual dye and ethidium bromide. This kit also allows for concentration of DNA, changing of buffers and desalting. Buffer PCR contains a pH indicator for easy determination of the optimal pH for efficient recovery of DNA fragments but at the same time does not interfere with DNA yield and quality as it is thoroughly removed during the subsequent washing steps. High recovery of pure DNA obtained is ready-to-use in many routine molecular biology applications such as restriction enzyme digestion, radioactive/fluorescence DNA sequencing, PCR, ligation and transformation, probe preparations and other manipulations.

### Features:

- Indicator dye for easy pH determination
- 90% of recovery achievable
- Mini-column spin technology
- Purification process less than 15 minutes

### Application:

For purification of DNA from reaction solutions

### Kit Components:

- GF-1 columns
- Collection tubes
- PCR DNA Binding Buffer (Buffer PCR)
- Wash Buffer (concentrate)\*
- Elution Buffer
- Handbook
- \* Please refer to Reconstitution of Solutions and Storage and Stability before using this kit.

Cat.-no	Description	Amount
PC05	PCR clean up extraction kit	5 preps
PC050	PCR clean up extraction kit	50 preps
PC100	PCR clean up extraction kit	100 preps
PC200	PCR clean up extraction kit	200 preps

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# Gel Extraction Kit

## Description:

The GF-1 Gel extraction Kit is a system designed for rapid purification of DNA bands ranging from 100bp to 10kb from all grades of agarose gel in TAE (Tris-acetate/ EDTA) or TBE (Tris-borate/ EDTA) buffer. Special buffers provided in the kit are optimized to enhance binding of DNA onto a specially-treated glass filter membrane for efficient recovery of highly pure DNA.

High recovery of pure DNA is obtained and ready to use in many routine molecular biology applications such as restriction enzyme digestion, radioactive/ fluorescence DNA sequencing, PCR, ligation, probe preparations and other manipulations

## Features:

- 90% of recovery achievable
- Mini-column spin technology
- Purification process less than 20 minutes
- Highly pure DNA ready to use for routine molecular biology applications such as restriction enzyme digestion, PCR, ligation, DNA sequencing, probe preparations, etc.

## Application:

Recovery/isolation of DNA from agarose gel

## Kit Components:

- GF-1 columns
- Collection tubes
- Gel DNA Binding Buffer (Buffer GB)
- Wash Buffer (concentrate)\*
- Elution Buffer
- Handbook

\* Please refer to Reconstitution of Solutions and Storage and Stability before using this kit.

Cat.-no	Description	Amount
GP05	Gel extraction kit	5 preps
GP050	Gel extraction kit	50 preps
GP100	Gel extraction kit	100 preps

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## AmbiClean Gel/PCR Purification

### Description:

The GF-1 AmbiClean Kit (Gel & PCR) contains special buffers to provide the correct salt concentration and pH for efficient recovery (80 - 90%) of DNA from both PCR product and agarose gel from both TAE or TBE buffer. This "clean up" - kit uses a specially treated glass filter membrane fixed into a column to efficiently bind DNA in the presence of high salt. It applies the principle of a mini-column spin technology and is well suited for the removal of agarose, excess dNTPs, short oligo fragments, mineral oil, enzymes from a PCR reaction product, proteins after restriction enzyme treatment and dephosphorylation and residual dye and ethidium bromide. This extraction kit also allows for concentration of DNA, changing of buffers and desalting. High recovery of pure DNA obtained is ready-to-use in many routine molecular biology applications such as restriction enzyme digestion, radioactive/fluorescence DNA sequencing, PCR, ligation and transformation, probe preparations and other manipulations.

### Kit Components:

- GF-1 columns
- Collection tubes
- DNA Binding Buffer (Buffer DB)
- Wash Buffer (concentrate)\*
- Elution Buffer
- Handbook

\* Please refer to Reconstitution of Solutions and Storage and Stability before using this kit.

Cat.-no	Description	Amount
GV05	AmbiClean PCR/Gel DNA extraction kit	5 preps
GV050	AmbiClean PCR/Gel DNA extraction kit	50 preps
GV100	AmbiClean PCR/Gel DNA extraction kit	100 preps

## RNA Purification Kits

- Total RNA Extraction Kit ..... 28-29
- Blood Total RNA Extraction Kit ..... 30-31
- Viral Nucleic Acid Extraction Kit ..... 32-33

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# Total RNA Extraction Kit

## Description:

The GF-1 Total RNA Purification Kit is designed for rapid and efficient purification of total RNA from various types of samples including cultured animal cells and tissues. The purification is based on the usage of denaturing agents to provide efficient cell lysis, denaturation of proteins and subsequent release of RNA. Special buffers provided in the kit are optimized to enhance the binding of RNA onto a specially treated glass filter membrane for efficient recovery of highly pure RNA.

## Features:

- Yields up to 100 µg of total RNA.
- Mini-column spin technology.
- No organic-based extraction required.
- Highly pure total RNA, ready to use for RT-PCR, Northern Blotting, polyA RNA (mRNA) purification, nuclease protection and in vitro translation.

## Application:

For purification of total RNA from cell based samples

## Kit Components:

- RNA binding columns
- Homogenization columns
- Collection tubes
- Buffer TR\*
- Inhibitor Removal Buffer\*
- Wash Buffer\*
- DNase I\*
- Digestion Buffer
- Digestion Enhancer
- RNase-free Water
- Proteinase K\*
- Handbook

\* Please refer to Reconstitution of Solutions and Storage and Stability before using this kit.

Cat.-no	Description	Amount
TR05	Total RNA Extraction Kit	5 preps
TR050	Total RNA Extraction Kit	50 preps
TR100	Total RNA Extraction Kit	100 preps

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## Total RNA Blood extraction

### Description

The GF-1 Total Blood RNA Extraction Kit is designed for rapid and efficient purification of total RNA from up to 1 ml fresh or frozen anti-coagulated whole blood. The purification is based on the usage of denaturing agents to provide efficient cell lysis, denaturation of proteins and subsequent release of RNA. Special buffers provided in the kit are optimized to enhance the binding of RNA onto a specially-treated glass filter membrane for efficient and fast recovery of highly pure RNA.

### Features

- Mini-column spin technology.
- Yields up to 4µg of total RNA.
- No organic-based extraction required.
- Highly pure total RNA, ready to use for RT-PCR, Northern Blotting, polyA RNA (mRNA) purification, nuclease protection and in vitro translation.

### Application

For purification of human total blood RNA

### Kit Components:

- Homogenization Columns
- RNA Binding Columns
- Collection tubes
- Buffer BR\*
- Inhibitor Removal Buffer\*
- Wash Buffer\*
- Proteinase K\*
- DNase I\*
- Digestion Buffer
- Digestion Enhancer
- RNase-free Water
- Handbook \* Please refer to Reconstitution of Solutions and Storage and Stability before using this kit.

Cat.-no	Description	Amount
TB05	Total RNA Blood extraction kit	5 preps
TB025	Total RNA Blood extraction kit	25 preps

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# Viral Nucleic Acid Extraction Kit

## Description:

The GF-1 Viral Nucleic Acid Extraction Kit is designed for rapid and efficient purification of viral DNA/RNA from samples such as serum, plasma, body fluid or virus-infected cell culture supernatant. The purification is based on the usage of denaturing agents to provide efficient viral lysis, denaturation of proteins and subsequent release of DNA or RNA. Special buffers provided in the kit are optimized to enhance the binding of DNA or RNA onto a specially-treated glass filter membrane for efficient recovery of highly pure DNA or RNA.

## Kit Components:

- GF-1 columns
- Collection tubes
- Buffer VL
- Wash Buffer 1 (concentrate)\*
- Wash Buffer 2 (concentrate)\*
- Carrier RNA\*
- Elution Buffer
- Handbook

\* Please refer to Reconstitution of Solutions and Storage and Stability.

Cat.-no	Description	Amount
RD05	Viral Nucleic Acid Extraction Kit	5 preps
RD050	Viral Nucleic Acid Extraction Kit	50 preps

● GeneON ●