

Oligo (dt)₁₅

Description:

- The primer is designed to initiate synthesis of a cDNA from total RNA in a reverse transcription reaction.
- For use in generating labeled cDNA for screening microarrays.
- The primer hybridizes to the poly(A) tail of mRNA.

Amount supplied: 30 µg (MW: 4501 mol/l; **Tm:** 26°C) = 1 OD unit (A260)

Storage: store at - 20°C

Transport: with "blue ice"

General Protocol:

Place fresh microtube in ice and make reaction mixture:

- **RNA template** 2 microliters
- Total RNA 0.1 - 5 micrograms
- or poly(A)+RNA 10 - 0.5 nanograms
- or specific RNA 0.01 picograms
- **random primer** 1 microliter (0.5 microgram)
- **RNAse-free water to** 16 microliters

Mix, collect by brief centrifugation.

2. Incubate mixture for 5 min at +70 °C, place in ice, collect by brief centrifugation.

3. Place microtube in ice and add:

- **x5 RT-buffer** 5 microliters
- **2.5 mM dNTP mixture** 4 microliters
- **M-MLV RT** 0.5 microliters

4. Incubate mixture for 10 min at +25 °C and 60 min at +37 °C .

5. Stop reaction by heating for min at +70 °C.

Ordering information:

Cat.-no	Description	Amount
S140	Oligo (dt)15	30 µg

GeneON .. a good decision ..

Contact Phone +49-(0)-621- 5720 864 Fax: +49-(0)-621-5724 462

E-Mail: <mailto:info@geneon.net> WEB: <http://www.taq-dna.com/> Version: 25.10.2009 AS

Unless specified otherwise, all products of GeneON are sold for research use only.

.. a good decision ..

Oligo (dt)₁₅
Cat: S140, 30µg

Description:

- The primer is designed to initiate synthesis of a cDNA from total RNA in a reverse transcription reaction.
- For use in generating labeled cDNA for screening microarrays.
- The primer hybridizes to the poly(A) tail of mRNA.

Amount supplied: 30 µg (MW: 4501 mol/l; **Tm:** 26°C) = 1 OD unit (A260)

Storage: store at - 20°C

Transport: with "blue ice"

General Protocol:

Place fresh microtube in ice and make reaction mixture:

- **RNA template** 2 microliters
- Total RNA 0.1 - 5 micrograms
- or poly(A)+RNA 10 - 0.5 nanograms
- or specific RNA 0.01 picograms
- **random primer** 1 microliter (0.5 microgram)
- **RNAse-free water to** 16 microliters

Mix, collect by brief centrifugation.

2. Incubate mixture for 5 min at +70 °C, place in ice, collect by brief centrifugation.

3. Place microtube in ice and add:

- **x5 RT-buffer** 5 microliters
- **2.5 mM dNTP mixture** 4 microliters
- **M-MLV RT** 0.5 microliters

4. Incubate mixture for 10 min at +25 °C and 60 min at +37 °C .

5. Stop reaction by heating for min at +70 °C.

Ordering information:

Cat.-no	Description	Amount
S140	Oligo (dt)15	30 µg

GeneON .. a good decision ..

Contact Phone +49-(0)-621- 5720 864 Fax: +49-(0)-621-5724 462

E-Mail: mailto:info@geneon.net WEB: <http://www.taq-dna.com/> Version: 25.10.2009 AS

Unless specified otherwise, all products of GeneON are sold for research use only.

.. a good decision ..